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Technology for Vacuum Systems

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Safety information!

CAUTION

- Ensure that repair is done only by a suitable trained and supervised technician. Obey local and national safety requirements. Ensure that the technician is familiar with the safety procedures which relate to the substances processed by the pumping system. Ensure that the device is decontaminated before repair and that you take adequate precautions to protect people from the effects of dangerous substances if contamination has occurred.

NOTICE

Read the **instructions for use** for the equipment carefully, especially sections "Safety informations" and "Use and operation". If necessary order "Instructions for use" from VACUUBRAND.

WARNING

If hazardous or corrosive substances have been used:

- ☞ The device will be contaminated with the process chemicals which have been pumped during operation. Ensure that the device is decontaminated before maintenance.
- ☞ Take adequate precautions to protect people from the effects of dangerous substances. Wear appropriate safety-clothing, do not inhale and avoid contact with skin.
- ☞ Before starting repair, isolate the device and other components from the vacuum system and the **electrical supply** so that they **cannot be operated accidentally**. Allow the device to cool, so that it is at safe temperature for maintenance work.
- ☞ Before starting maintenance, **wait two minutes** after isolating the equipment from mains to allow the capacitors to discharge.

NOTICE

Before starting repair check that the required parts are available and of the correct type before you start your work.

Check the operating sequence mentally on feasibility, safety requirements and consequences on safety and function of the equipment.

WARNING

- ☞ Use only genuine spare parts and accessories.
If using components of other manufacturers the function and the safety of the equipment may be restricted.
- ☞ Do not reuse damaged parts.

CAUTION

- **Check operability and safety after repair.**
In case of problems return the device to the factory for inspection or repair if necessary.
- Obey regulations when disposing of solvents and chemicals.

NOTICE

In order to comply with law (occupational, health and safety regulations, safety at work law and regulations for environmental protection) vacuum pumps, components and measuring instruments returned to the manufacturer can be repaired only when certain procedures (see section "Notes on return to the factory" in the "Instructions for use") are followed.

Scrapping and waste disposal

Dispose of the pump and any components removed from it safely in accordance with all local and national safety and environmental requirements.

Particular care must be taken with components and waste oil which have been contaminated with dangerous substances from the process.

- ☞ You may authorize us to dispose the equipment at your expense.

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Upgrade kit 20699906:**Fitting the inlet of NT vacuum systems with a manometer at the inlet.**

The manometer is to be mounted at the distribution head of NT vacuum systems. The manometer may be mounted at the following vacuum systems:

MZ 2C NT + 2 AK
 MZ 2C NT + AK + EK
 MD 4C + 2 AK
 MD 4C + AK + EK

➔ **Before starting assembly work switch off the diaphragm pump and separate it from mains! Remove catchpot.**

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



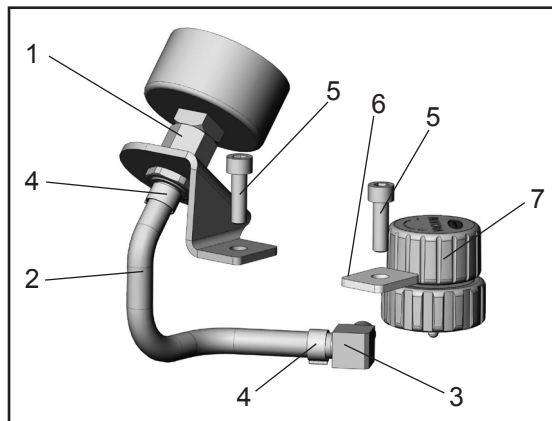
- 1: Manometer with mounting angle
- 2: Preformed hose
- 3: Connection angle
- 4: Hose clip
- 5: Screw 25mm
- 6: Height compensation
- 7: Flow control valve

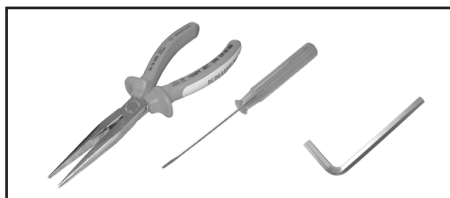Note: The upgrade kit contains surplus hose clips.



**Tools required:**

- hex key size 5
- flat pliers
- flat-bladed screw driver 2.5 mm

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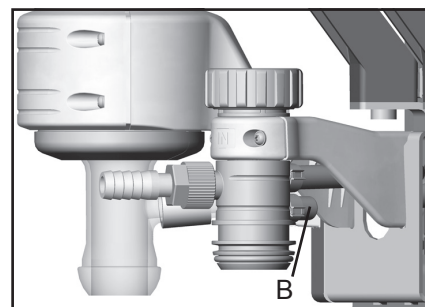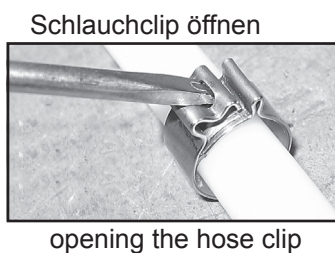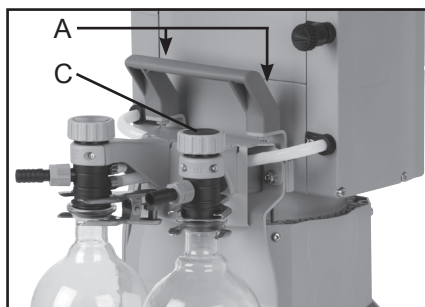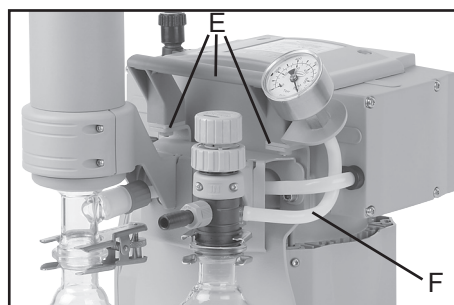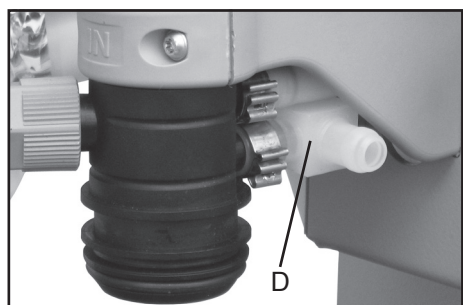Upgrade kit 20699907:

Fitting the inlet of NT diaphragm pumps/pumping units with a manometer at the inlet.

The manometer is to be mounted at the inlet block of NT pumps/pumping units. The manometer may be mounted at the following pumps/pumping units:

MZ 2C NT + AK SYNCHRO + EK
 MD 4C NT + AK SYNCHRO + EK
 PC 511 NT (at inlet with manual flow control valve)
 PC 611 NT (at inlet with manual flow control valve)

➔ **Before starting assembly work switch off the diaphragm pump and separate it from mains!**

- ➔ Unscrew the hose nozzle DN 10 mm at the inlet block. If necessary, use universal pliers carefully to loosen the hose nozzle.
- ➔ Screw the new hose nozzle with manometer into the thread at the inlet block. Pay attention to the seal ring.
- ➔ Align the manometer and tighten the assembly with a hex key size 6 only hand-tight. Attention: Do not overwind! (Thread is cut into plastics.)

Inlet: hose nozzle DN 10mm;
 Wetted materials: stainless steel

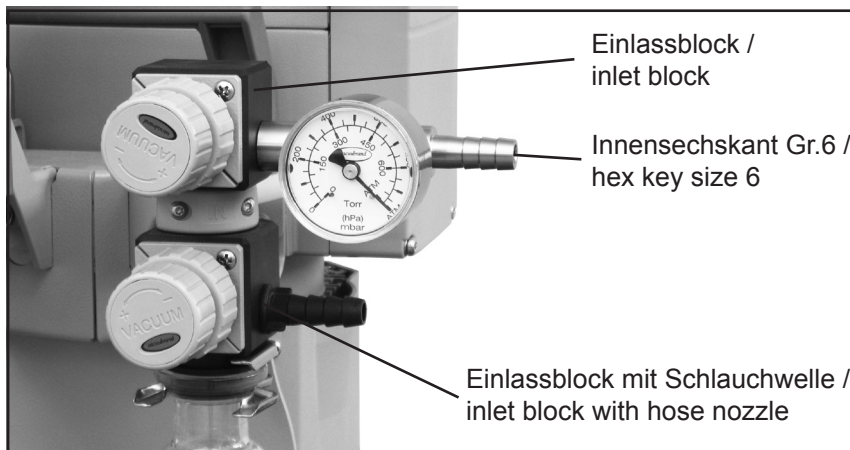
Abb.: Ausbausatz Manometer montiert an MD 4C NT +AK SYNCHRO + EK
 fig.: Upgrade kit manometer mounted at MD 4C NT + AK SYNCHRO + EK

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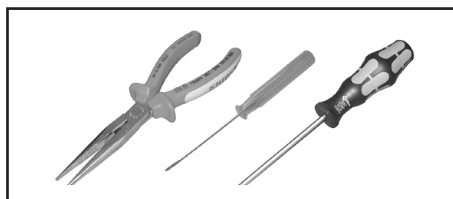
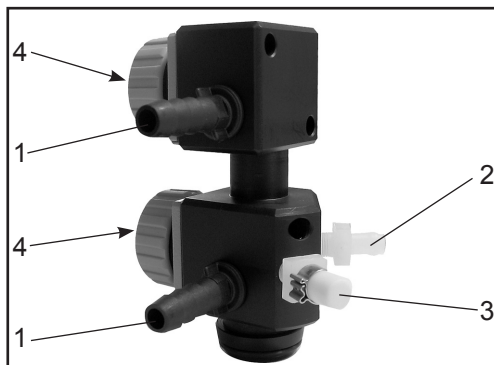
- 1: connection to application
 2: connection to pump
 3: measuring connection to controller
 4: blind plug

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Upgrade kit 2069920:**Fitting NT vacuum systems and pumping units mounted on metal support with two inlets.**

The upgrade kit is to be mounted at the inlet of the vacuum systems/pumping units. The upgrade kit may be mounted at the following vacuum systems/pumping units:

PC 510 NT
 PC 610 NT
 MZ 2C NT + 2 AK
 MZ 2C NT + AK + EK
 MD 4C NT + 2 AK
 MD 4C NT + AK + EK

➔ **Before starting assembly work switch off the diaphragm pump and separate it from mains! Remove catchpot.**

Required tools:

- Torx screw driver TX10
- flat-bladed screwdriver 2.5 mm
- flat pliers

Assembly work is shown on a pumping unit PC 510 NT. Carry out assembly work on other NT vacuum systems or pumping units in the same way. In case of vacuum systems without controller the measuring connection will remain unconnected.

1. Open one (vacuum systems) or two (pumping units) hose clips at the hoses at the inlet block (A): Open the hose clip with a flat-bladed screwdriver. Apply screw driver as shown in fig. B and turn.

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5. Assembling the new inlet block:
Only pumping units with controller:
Shorten the hose to the controller by 15mm. Put a hose clip onto the connection hose and connect the hose at the hose connection next to the hose nozzle.

Shorten the hose to the pump by 15mm. Put a hose clip onto the connection hose and push the hose onto the corresponding hose connection of the new inlet block.

Close the hose clips with flat pliers.

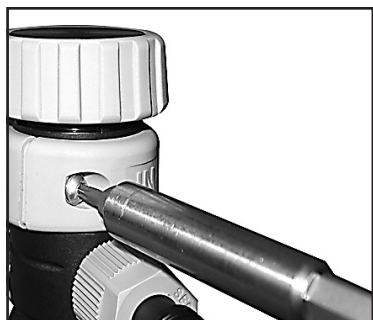
Assemble the inlet block with the counterholder at the the pumping unit (two Torx screws).

6. If using a solenoid operated in-line valve, remove the blind cap at the inlet block and assemble in-line valve.

Note: A solenoid operated in-line valve should be assembled at the lower inlet block (with measuring connection).

There is no measuring connection at the upper inlet block. If a solenoid valve at the upper inlet block shall be used, it is necessary to install a separate measuring connection to the controller, e. g. using a T-piece in the suction line.

Attention: The flow control valve at the upper inlet block and the solenoid operated valve at the lower inlet block shown in figure D are no included in the upgrade kit.

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Upgrade kit 20699942:**Fitting NT pumping units in plastic housing with two inlets.**

The upgrade kit is to be mounted at the inlet of the pumping units. The upgrade kit may be mounted at the following pumping units:

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
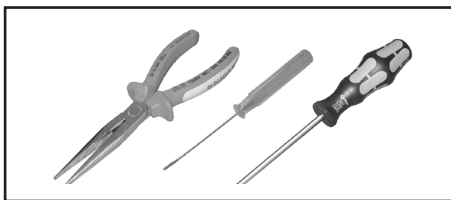
Assembly work is shown on a pumping unit PC 510 NT. Carry out assembly work on other NT pumping units in the same way.

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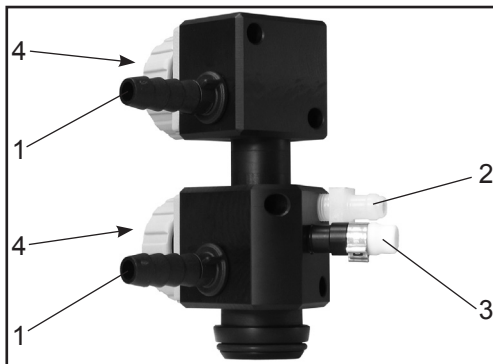
- 1: connection to application
- 2: connection to pump
- 3: measuring connection to controller
- 4: blind plug

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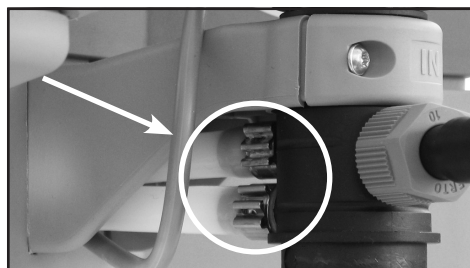1. Open the two hose clips at the hoses at the inlet block (A): Open the hose clip with a flat-bladed screwdriver. Apply screw driver as shown and turn.

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© 2018 VACUUBRAND GMBH + CO KG Printed in Germany