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VACUU·PURE 10
VACUU·PURE 10C

Montageanleitung Mounting instructions

Original instructions Keep for further use!

This manual is only to be used and distributed in its complete and original form. It is strictly the user's responsibility to carefully check the validity of this manual with respect to the product.

Manufacturer:

VACUUBRAND GMBH + CO KG
Alfred-Zippe-Str. 4
97877 Wertheim
GERMANY

Phone:

Head office +49 9342 808-0
Sales +49 9342 808-5550
Service +49 9342 808-5660

Fax: +49 9342 808-5555

Email: info@vacuubrand.com

Web: www.vacuubrand.com

*Thank you for purchasing this product from **VACUUBRAND GMBH + CO KG**. You have chosen a modern and technically high quality product.*

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6 Introduction

This installation manual is part of your product. The installation manual describes how to replace spare parts for the vacuum pumps *VACUU-PURE 10* and *VACUU-PURE 10C*. The installation manual applies to all versions of the pump and is intended in particular for specialists.

6.1 User information

Safety

Installation manual
and safety

- Read this installation manual thoroughly and completely before starting maintenance work.
- In addition to the information in this manual, adhere to the accident prevention regulations and industrial safety regulations applicable in the country of use.

General

General
information

- To improve readability, the general term *vacuum pump* is used as an equivalent to and instead of the product name.
- The illustrations in this manual are only intended to facilitate comprehension.
- We reserve the right to make technical changes in the course of continuous product improvement.

Copyright

Copyright © and
copyright law

The content of this manual is protected by copyright. Only copies for internal use are allowed, e.g., for professional training.

© VACUUBRAND GMBH + CO KG

Contact

Contact
us

- If your manual is incomplete, you can request a replacement. Alternatively, you can use our download portal: www.vacuubrand.com
- You are welcome to contact us at any time in writing or by telephone if you would like more information, have questions about our products or wish to share feedback with us.
- When contacting our Service Department, please have the serial number and product type at hand → *see the rating plate on the product.*

6.2 About this document

6.2.1 Display conventions

Warning levels

Display conventions

	WARNING
	<p>Warns of a potentially hazardous situation. Disregarding the situation could result in serious injury or death.</p> <p>⇒ Take appropriate action to avoid dangerous situations!</p>

	CAUTION
	<p>Indicates a potentially hazardous situation. Disregarding the situation could result in minor injury or damage to property.</p> <p>⇒ Take appropriate action to avoid dangerous situations!</p>

NOTE
<p>Indicates a potentially harmful situation. Disregarding the situation could result in damage to property.</p>

Additional notes

IMPORTANT!

- ⇒ Information or specific recommendation which must be observed.
- ⇒ Important information for trouble-free operation of your product.

6.2.2 Symbols and icons

This manual uses symbols and icons. Safety symbols indicate specific risks associated with handling the product. Symbols and icons are designed to help you identify risks more easily.

Safety symbols

Explanation of safety symbols

Hazardous substance – hazards to human health.

Danger: electricity.

Warning: hot surface.

Disconnect power plug.

Wear chemical-resistant protective gloves.

Wear protective goggles.

Additional symbols and icons

Additional symbols

Positive example – **Do this!**
Result – **OK**

Negative example – **Don't do this!**

Refers to content in other supplementary documents.

6.2.3 Handling instructions (action steps)

Instructions (single step)

Instructions

⇒ Perform the step described.

Result of action

Instructions (multiple steps)

1. First step

2. Next step

Result of action

Perform the steps in the order described.

7 Safety information

The information in this chapter must be observed by everyone who works with the product described here.

Usage Only use the product if it is in perfect and complete working condition.

- Observe the installation manual
- Observe the vacuum pump instructions for use
- Have the vacuum pump inspected regularly according to its operating conditions by qualified personnel
- Use only approved spare parts

In the case of activities which require protective clothing, personal protective equipment as specified by the operator is to be worn.

If the product is not in proper working order, it must be prevented from being accidentally switched back on.

- ⇒ Always be conscious of safety and work in a safe manner.
- ⇒ Observe instructions issued by the operator, and national regulations on accident prevention and industrial safety.

7.1 Target group description

The installation manual must be read and observed by every person who is tasked with the activities described below.

Personnel qualification

Operator	Laboratory staff, such as chemists, laboratory technicians
Specialist	Person with professional qualification in mechanics, electrical equipment or laboratory devices
Responsible specialist	Specialist with additional specialist, departmental or area responsibility

Responsibility matrix

Activity	Operator	Specialist	Responsible specialist
Replace filters		x	x
Replace O-ring	x	x	x
Replace flat gasket	x	x	x
Replace blind plug	x	x	x
Shutdown		x	x
Repair order			x

7.2 Safety

IMPORTANT!

- ⇒ Only use original spare parts. The validity of the CE mark may be invalidated if original parts are not used.
- ⇒ When handling contaminated parts, follow the relevant regulations and safety precautions.

7.3 Protective clothing

When performing maintenance work, we recommend wearing chemical-resistant protective gloves, protective clothing and protective goggles, if there is a possibility that you could come into contact with contaminated parts.

IMPORTANT!



- ⇒ When handling chemicals, wear your personal protective equipment.

8 Spare parts

The following spare parts are available for the vacuum pump:

Overview of spare parts

Overview of spare parts

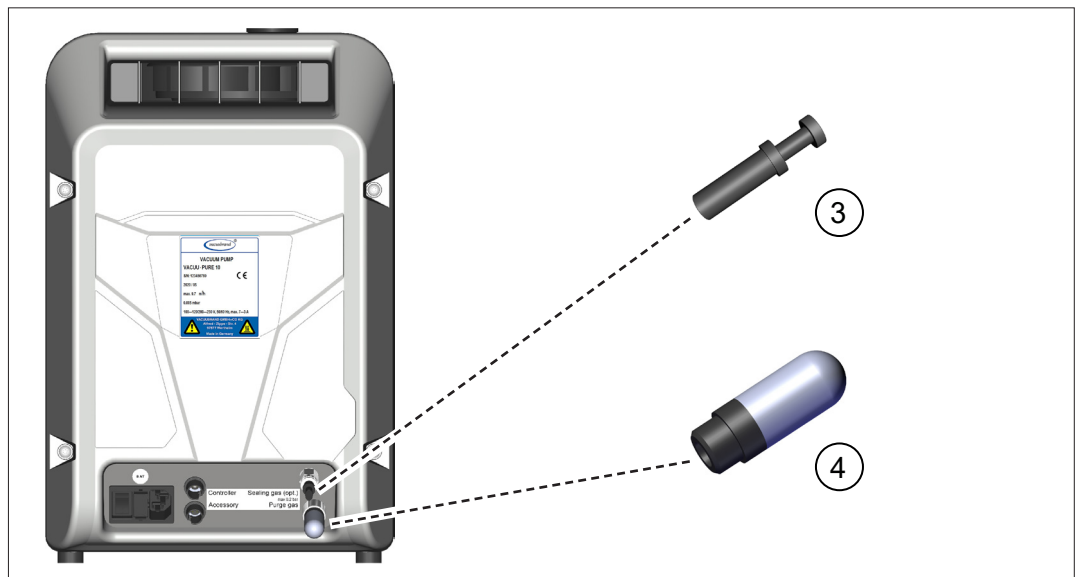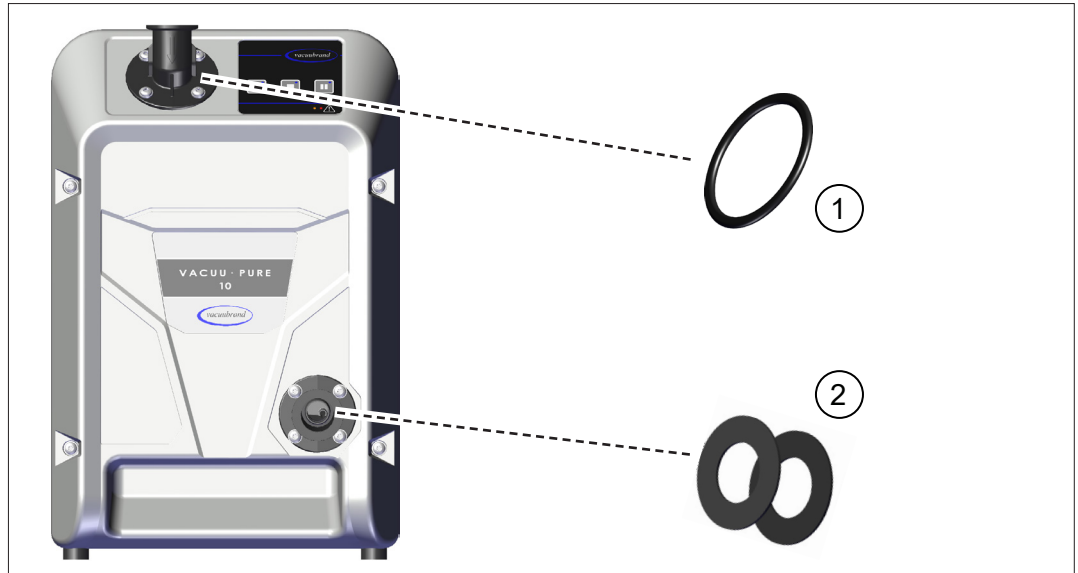Description

1	O-ring on the inlet flange	20638419
2	Flat gasket at the outlet	20638420
3	Blind plug, only VACUU·PURE 10	20638414
4	Air inlet filter, VACUU·PURE 10C (plus another 2 in the pump interior)	20638411

9 Replace spare parts

Replace dirty or damaged parts.

The vacuum pump instructions for use describe how to replace the air inlet filter for regeneration mode and how to replace the device fuses.

	WARNING
	Danger due to electrical voltage. ⇒ Switch the device off before cleaning or maintenance work. ⇒ Unplug the power plug from the socket.

NOTE
Damage possible if work is performed incorrectly. ⇒ Have maintenance work performed by a trained professional or at least by a trained person. ⇒ Recommendation: Before carrying out maintenance work for the first time, please read through all the handling instructions to get an overview of the required service work.

Tool Required tools

- Screwdriver TX25

9.1 Air inlet filter (#20638411)


Air inlet filter for the internal sealing gas supply of the VACUU·PURE 10C

The air for the internal sealing gas supply of the VACUU·PURE 10C is aspirated through two filters. Replace dirty or blocked air inlet filters for the internal sealing gas supply.


	CAUTION
	<p>Risk of burns from hot surfaces. During operation, internal parts of the pump can heat up to temperatures > 65 °C.</p> <ul style="list-style-type: none"> ⇒ Allow the vacuum pump to cool down before starting maintenance work. ⇒ Allow the vacuum pump to cool down before removing housing sections.

Remove dirty air filters


Remove air filter at rear of device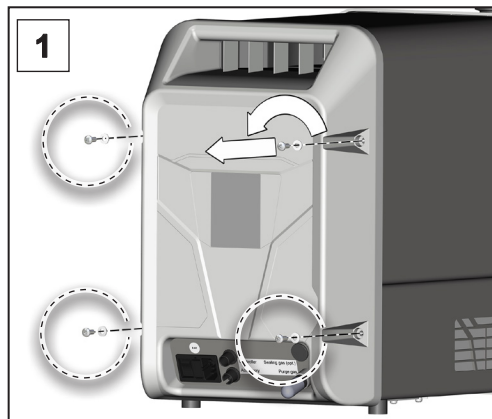


1. Undo the 4 screws on the rear of the housing using a Torx screwdriver TX25, paying attention to the washers.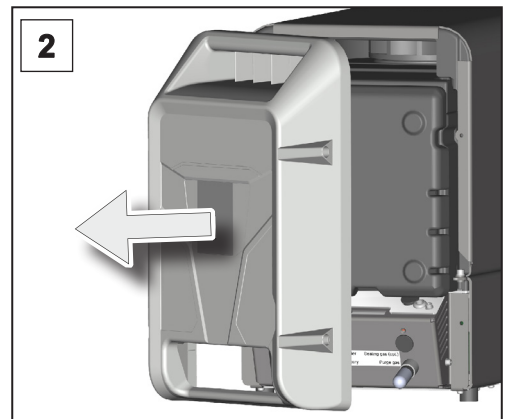

2. Remove the back of the housing.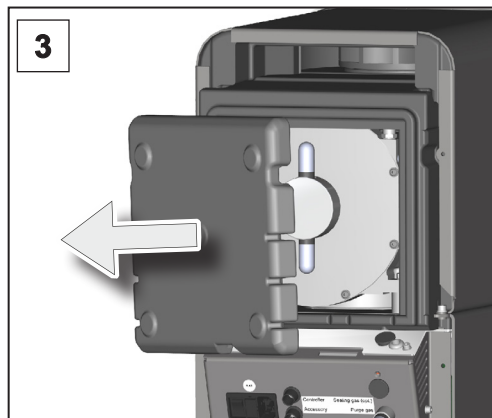


3. Remove the sound insulation.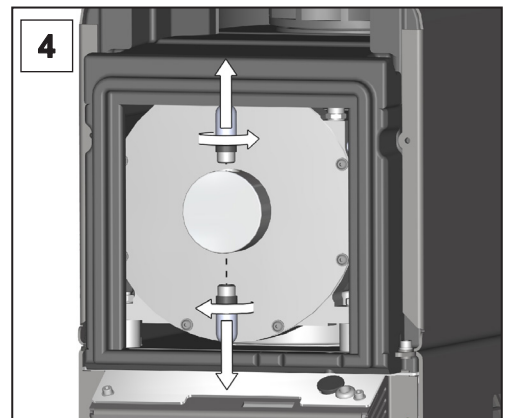


4. Unscrew the two dirty filters from the sealing gas supply inlet.

Install new air filters


Install air filters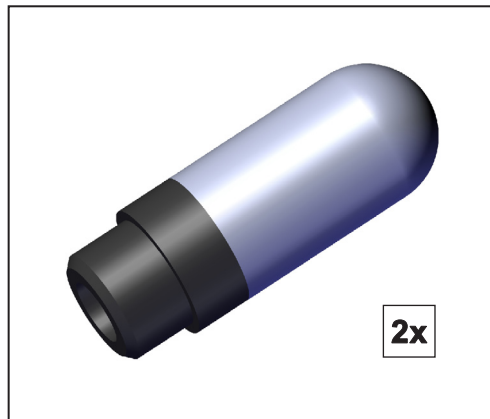


Filters (2x) in the inlet of the sealing gas supply.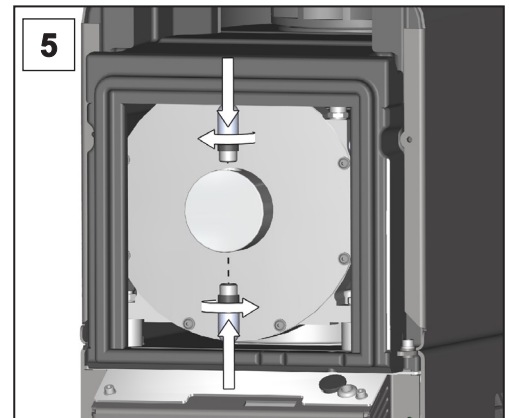


5. Screw the two new filters into the sealing gas supply inlet.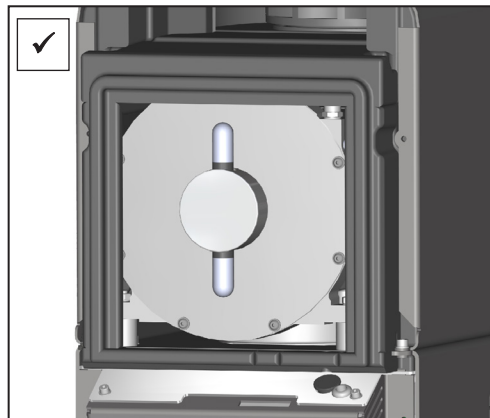


Both filters at the inlet of the sealing gas supply have been replaced.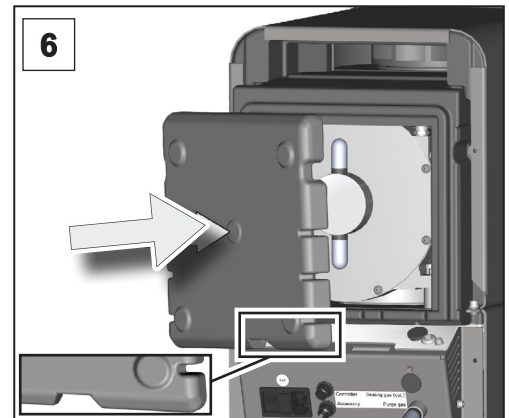

6. Refit the sound insulation on the back of the pump with the cable recess at the bottom.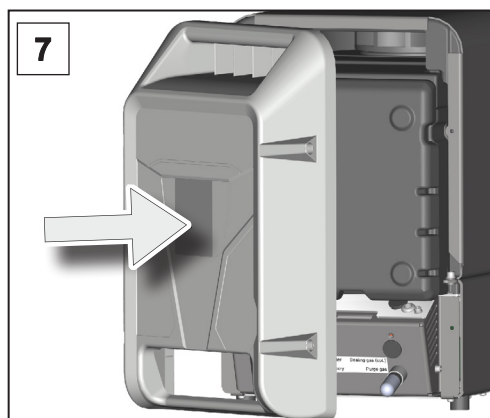

7. Refit the back of the housing.

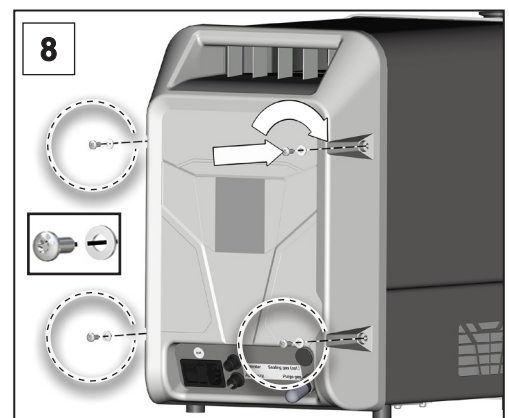8. Tighten the 4 screws on the rear of the housing until hand-tight using a Torx screwdriver TX25, paying attention to the washers.

9.2 O-ring on the inlet flange (#20638419)


Replace O-ring at inlet

Replace the O-ring on the inlet flange if dirty or damaged.


WARNING	
	<p>Risk from contaminated parts. Pumping hazardous media can result in hazardous substances adhering to internal parts of the pump.</p>
	<p>⇒ Wear your personal protective equipment, e.g., protective gloves, eye protection and, if necessary, respiratory protection.</p>
	<p>⇒ Decontaminate wetted parts that you may come into contact with during maintenance work.</p>
	<p>⇒ Take safety precautions according to the instructions you have received on handling hazardous substances.</p>

Remove the O-ring


Remove O-ring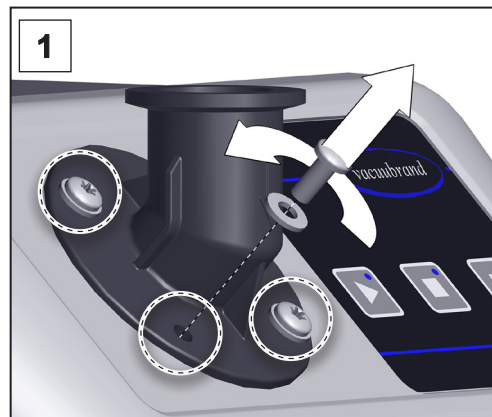

1. Undo the 4 screws on the inlet flange using a Torx screwdriver TX25, paying attention to the washers.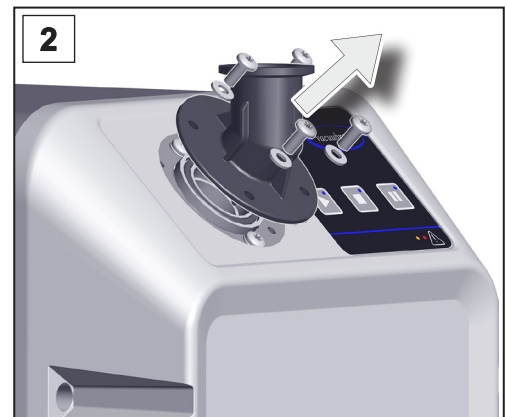


2. Remove the inlet flange.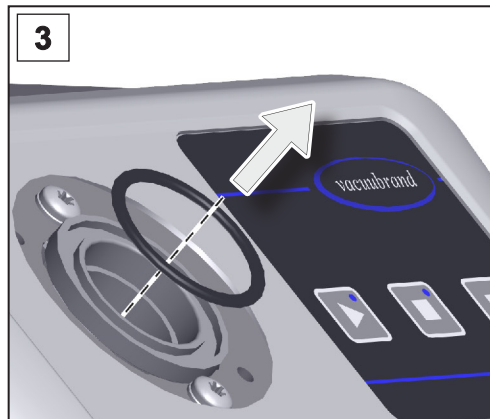

3. Remove the O-ring. Check the O-ring for damage and dirt.

Insert a new O-ring


Insert O-ring

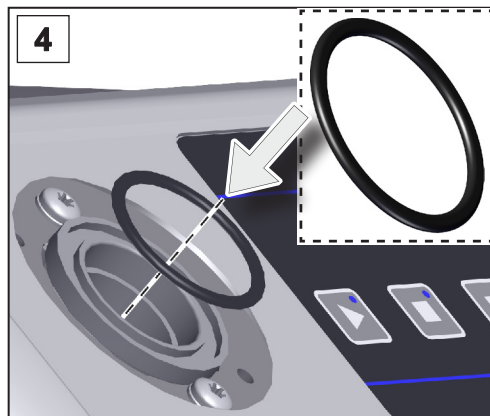4. Replace the O-ring. Insert the new O-ring.

5. Refit the inlet flange, paying attention to the washers: curved side towards the screw head.

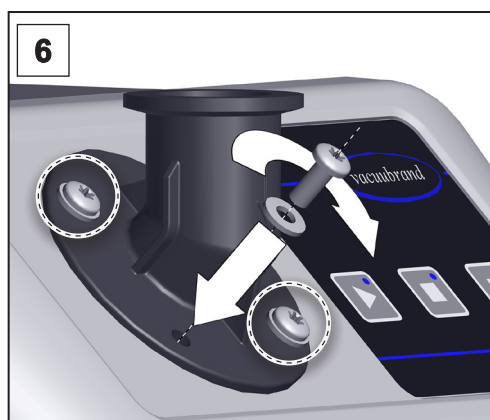6. Screw in the 4 screws on the inlet flange using a Torx screwdriver TX25.

9.3 Flat gasket at the outlet (#20638420)

Replace flat gasket at the outlet


Replace the flat gasket on the outlet connector if dirty or damaged.

	WARNING
 	<p>Risk from contaminated parts. Pumping hazardous media can result in hazardous substances adhering to internal parts of the pump.</p> <ul style="list-style-type: none"> ⇒ Wear your personal protective equipment, e.g., protective gloves, eye protection and, if necessary, respiratory protection. ⇒ Decontaminate wetted parts that you may come into contact with during maintenance work. ⇒ Take safety precautions according to the instructions you have received on handling hazardous substances.


	CAUTION
	<p>Risk of burns from hot surfaces. During operation, internal parts of the pump can heat up to temperatures > 65 °C.</p> <ul style="list-style-type: none"> ⇒ Allow the vacuum pump to cool down before starting maintenance work. ⇒ Allow the vacuum pump to cool down before removing housing sections.

Remove the flat gasket


Remove the flat gasket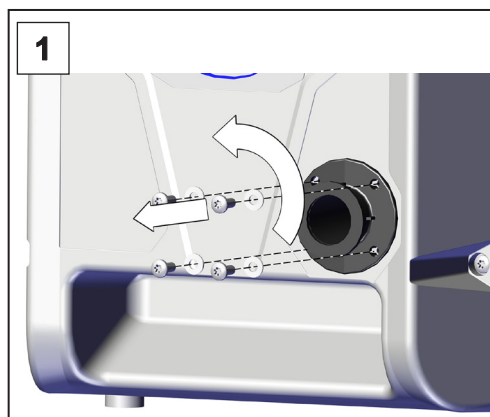


1. Undo the 4 screws on the outlet flange using a Torx screwdriver TX25, paying attention to the washers.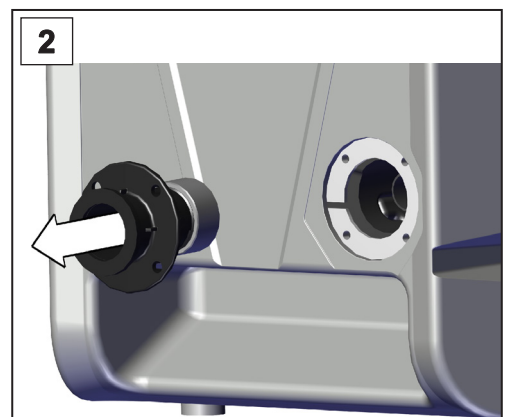

2. Remove the outlet flange.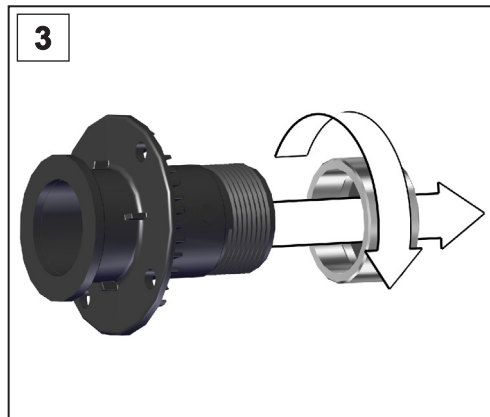


3. Unscrew the union nut from the outlet connector.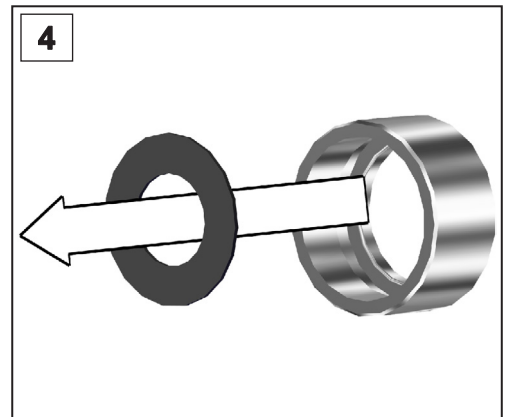


4. Remove the flat gasket. Check the gasket for damage and dirt.

Insert the flat gasket


Insert the flat gasket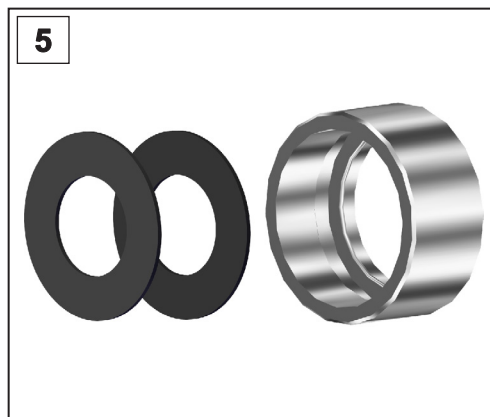


5. The gasket consists of two individual sealing membranes. Replace damaged or dirty gaskets.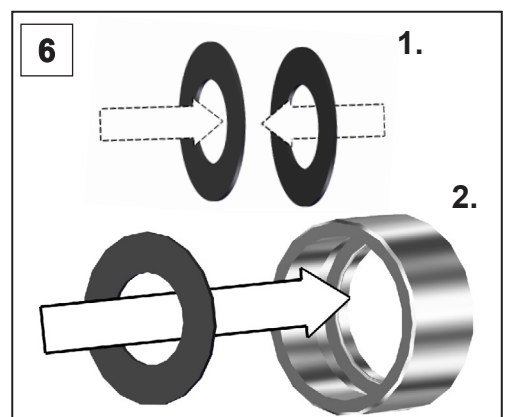


6. Put the two sealing membranes together (1.) and insert them into the union nut (2.).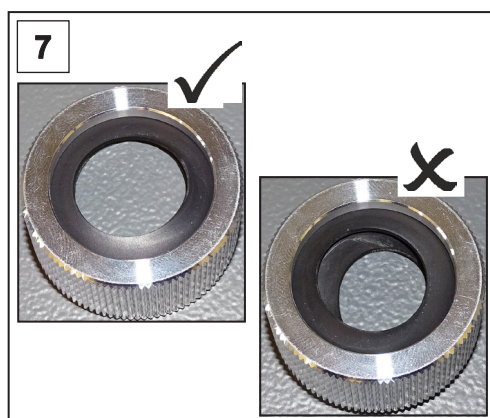


7. Ensure the sealing membranes are correctly positioned in the union nut.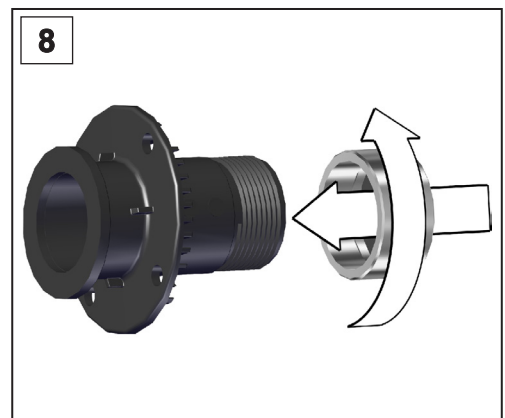


8. Screw the union nut with the gasket back onto the outlet connector until hand-tight.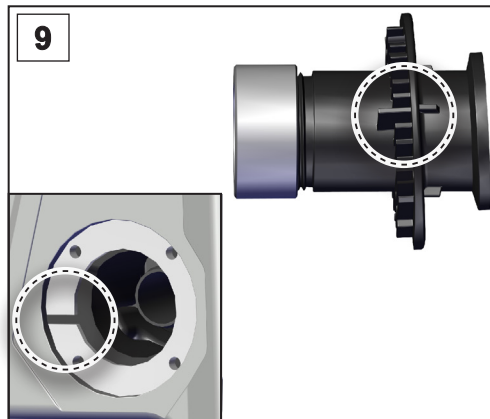

9. When positioning the outlet flange, ensure the protrusion on the outlet flange clicks into place in the recess of the pump.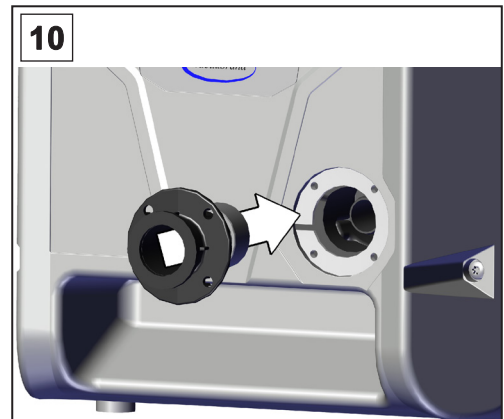

10. Refit the outlet connector.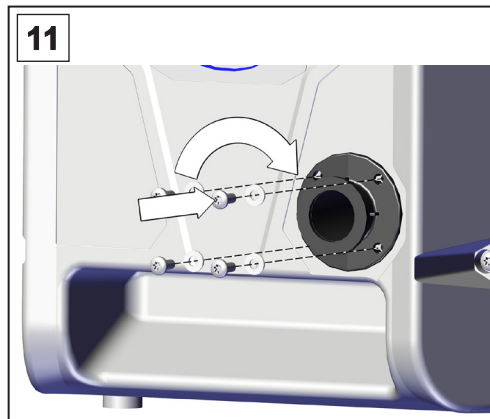

11. Screw in the 4 screws on the inlet flange using a Torx screwdriver TX25, paying attention to the washers: curved side towards the screw head.


9.4 Blind plug (#20638414)

Blind plug of
VACUU·PURE 10

Replace the blind plug on the VACUU·PURE 10 if damaged or missing.

Remove the blind plug


Remove the blind
plug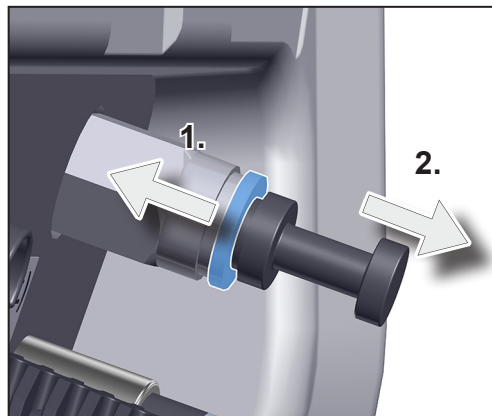

⇒ Press the blue ring towards the vacuum pump (1.) and take out the blind plug (2.).

Insert the blind plug

Insert the blind plug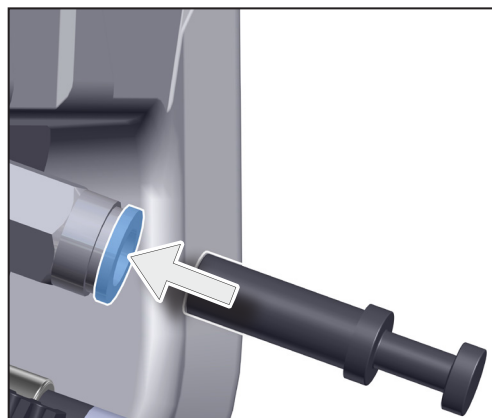

⇒ Push in the blind plug as far as it will go.

10 Appendix

Ordering information

Ordering information
for spare parts

Spare parts	Order no.
Air inlet filter (regeneration mode, internal sealing gas supply of the VACUU·PURE 10C)	20638411
Device fuses 8 AT 5x20	20638416
O-ring inlet flange	20638419
Flat gasket at outlet (FKM)	20638420
Blind plug (VACUU·PURE 10)	20638414


Sources of supply

Purchase original accessories and original spare parts from a subsidiary of **VACUUBRAND GMBH + CO KG** or your local distributor.

Information about our complete product range is available in the current [product catalog](#).

⇒ Your local distributor or **VACUUBRAND GMBH + CO KG** [sales office](#) is available to assist you with orders, questions on vacuum control and optimal accessories.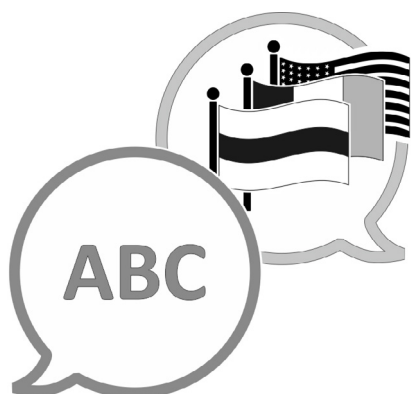

[VACUUBRAND > Support > Manuals](#)

Manufacturer:

VACUUBRAND GMBH + CO KG
Alfred-Zippe-Str. 4
97877 Wertheim
GERMANY

Phone:

Head office +49 9342 808-0
Sales +49 9342 808-5550
Service +49 9342 808-5660

Fax: +49 9342 808-5555

Email: info@vacuubrand.com

Web: www.vacuubrand.com

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Copyright © und